




K7003-72 PARTS LIST			
ITEM	QTY	PART NUMBER	DESCRIPTION
1	2	C7002-L	SHELF BRACKET, BOTTOM, LH FRONT, UTILITY
2	2	C7002-R	SHELF BRACKET, BOTTOM, RH FRONT, UTILITY
3	1	C7003-72 OR C7003-48	SHELF, BOTTOM, WELDED, UTILITY 72x24 OR 48 x 24
4	8	H236	SAE WASHER, #8
5	8	H183	SCREW, 8-32 THRD-CUT, 90087A194_

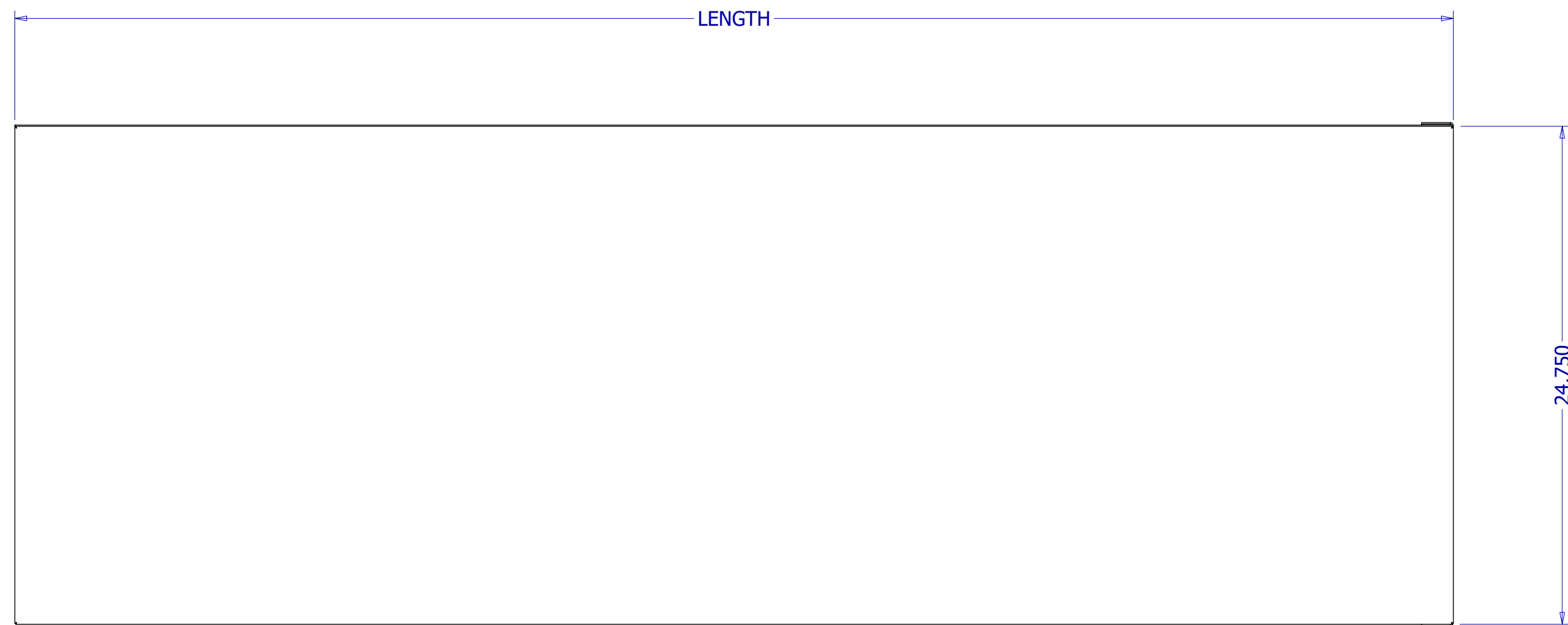

PART #	LENGTH
K7003-48	71.50"
K7003-72	59.50"

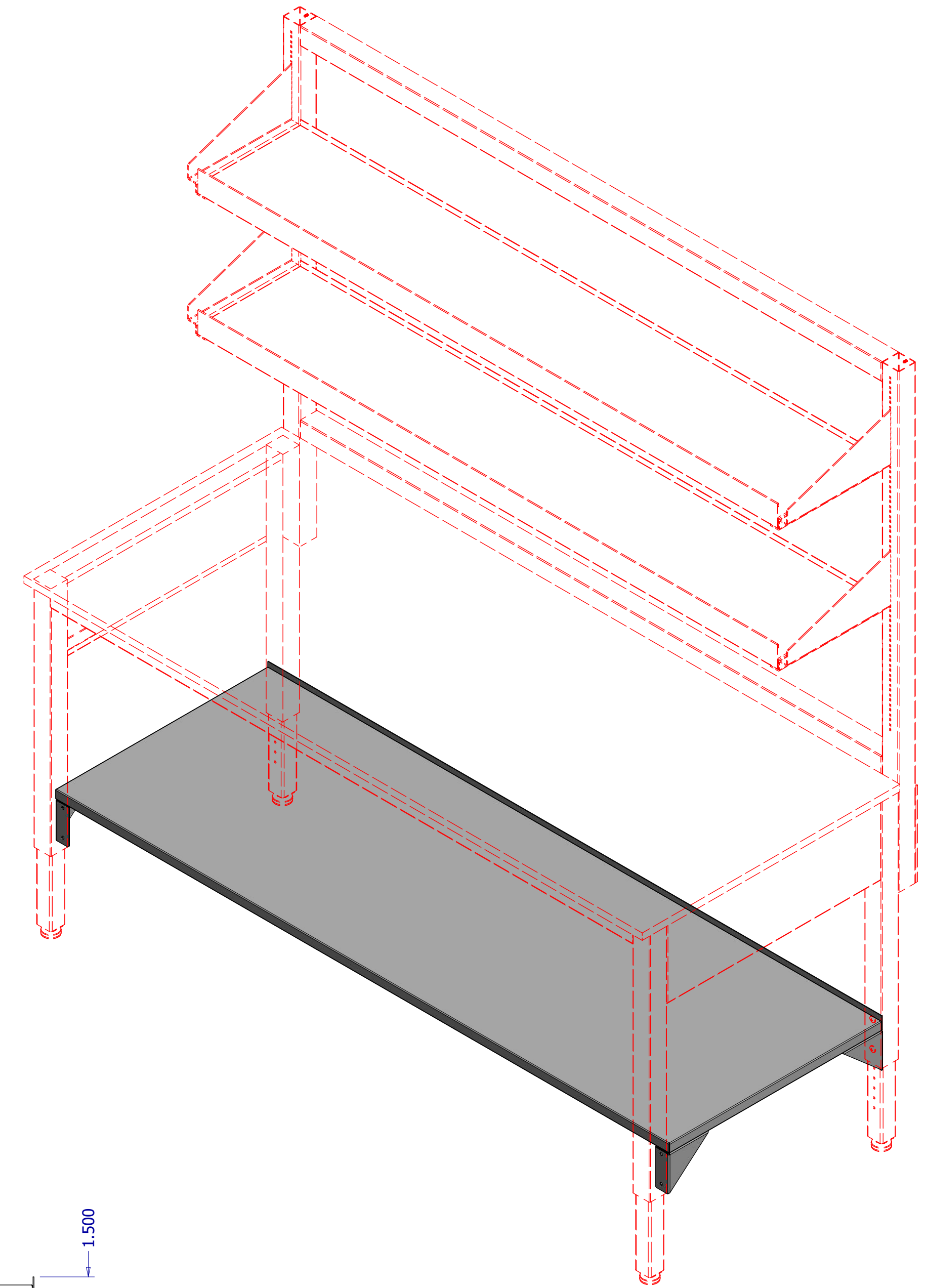

NOTE:
FIXED HEIGHT TABLES
REQUIRE MINOR DRILLING FOR
SHELF BRACKET ATTACHMENT

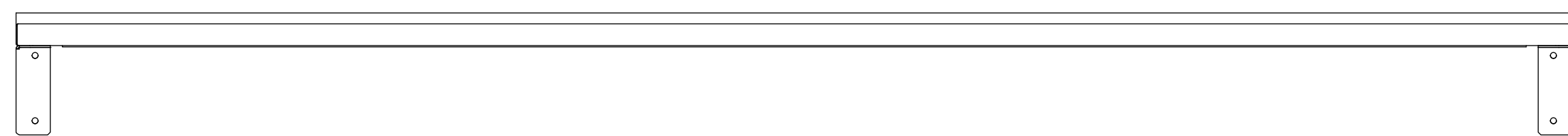

DRAWN AG	11/11/2020	LOOPED LOGIC 747 Calle Plano, Camarillo CA 93012 805-498-3121	
CHECKED		TITLE CUTSHEET LOOPED LOGIC LOWER SHELF KIT	
QA		FINISH PETAL WHITE	
APPROVED		SIZE D	PART NO
UNLESS OTHERWISE NOTED, GEN. TOLERANCE +/- .015"; ALL DIMENSIONS ARE IN INCHES	JOB # QUOTE #	SCALE 1/6" MATL	REV -
		SHEET 1 OF 1	