



BTUP48-G2


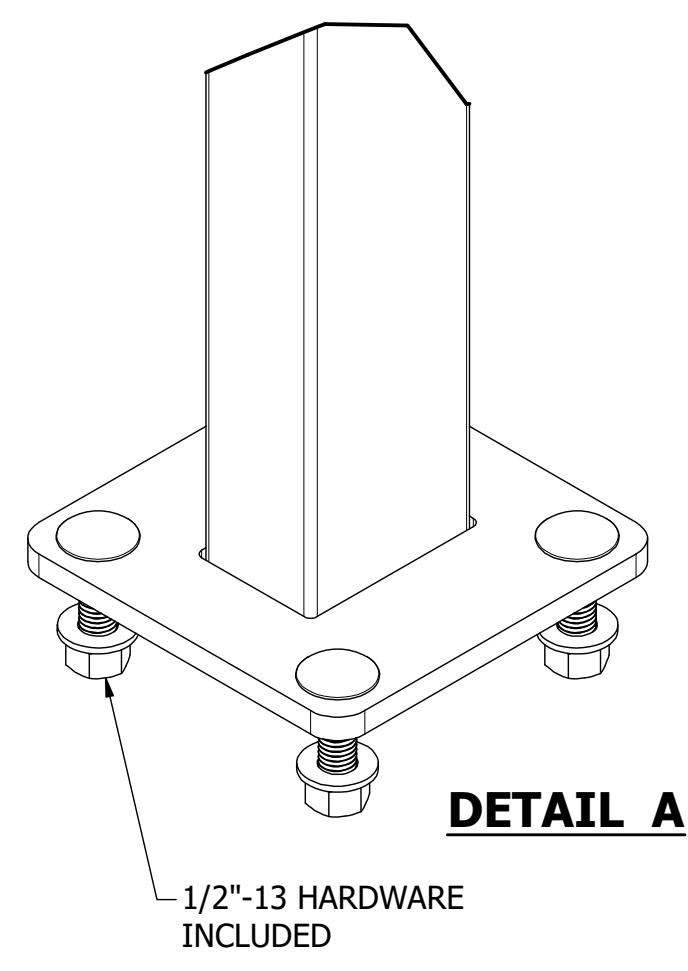


DETAIL A

1/2"-13 HARDWARE INCLUDED

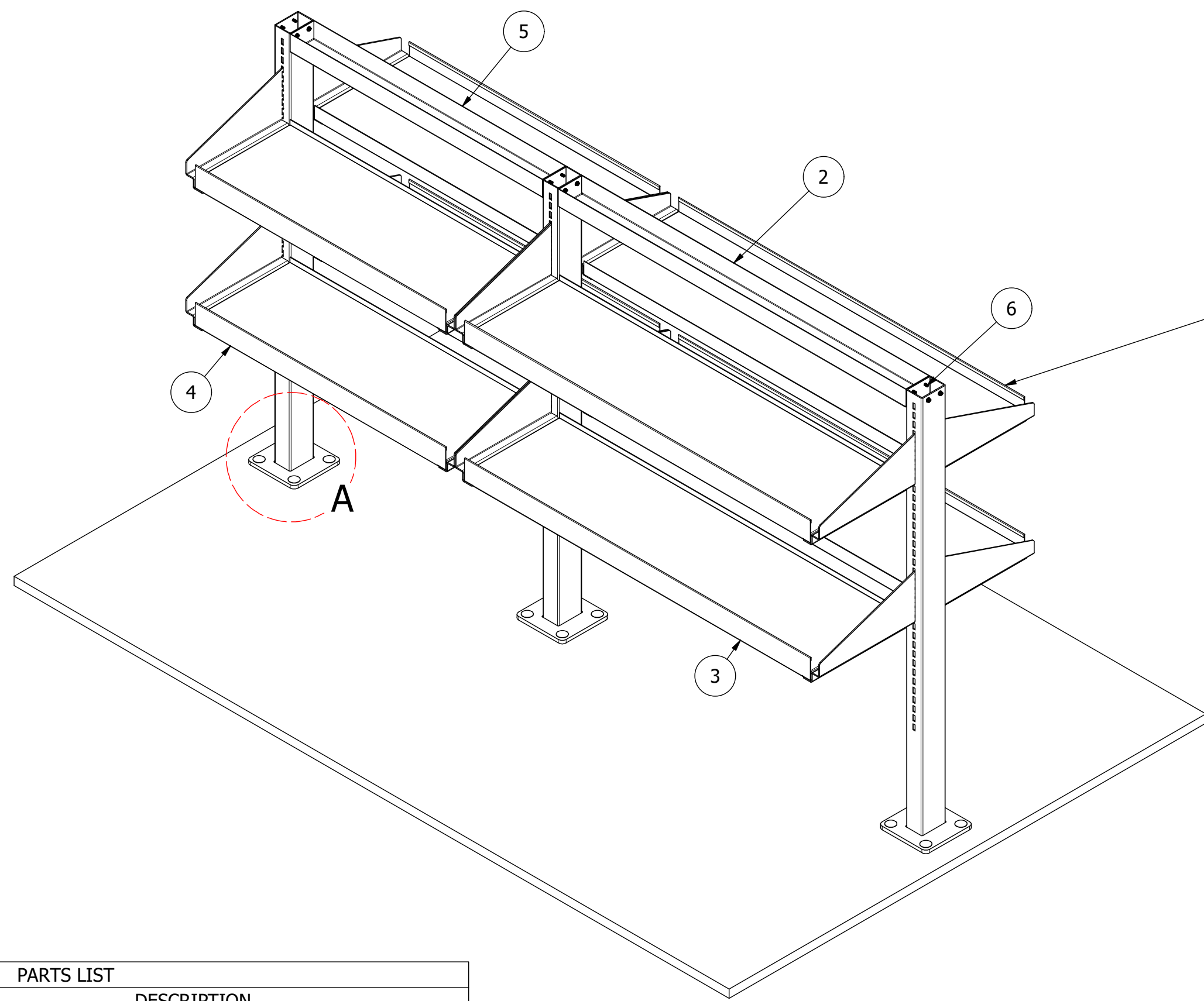

PARTS LIST			
ITEM	QTY	PART NUMBER	DESCRIPTION
1	3	BTUP48-G2	BENCH TOP MOUNTED UPRIGHT POST W/ MOUNT HARDWARE
2	1	C13174-43	HORIZONTAL UPRIGHT BRACE, FMUP, BTUP 2X3X43.75, G2
3	4	K13032-12	UPPER SHELF KIT WITH BRACKETS 48" x 12"
4	4	K13070-12	SHELF KIT, UPPER W BRKTS, 31.5X12, UTILITY TBL
5	1	C13174-31	BRACE, HORIZONTAL, FMUP/BTUP TOP RAIL, 2X3X31.75, G2
6	8	H237	SCREW, 10-24 THRD-CUT

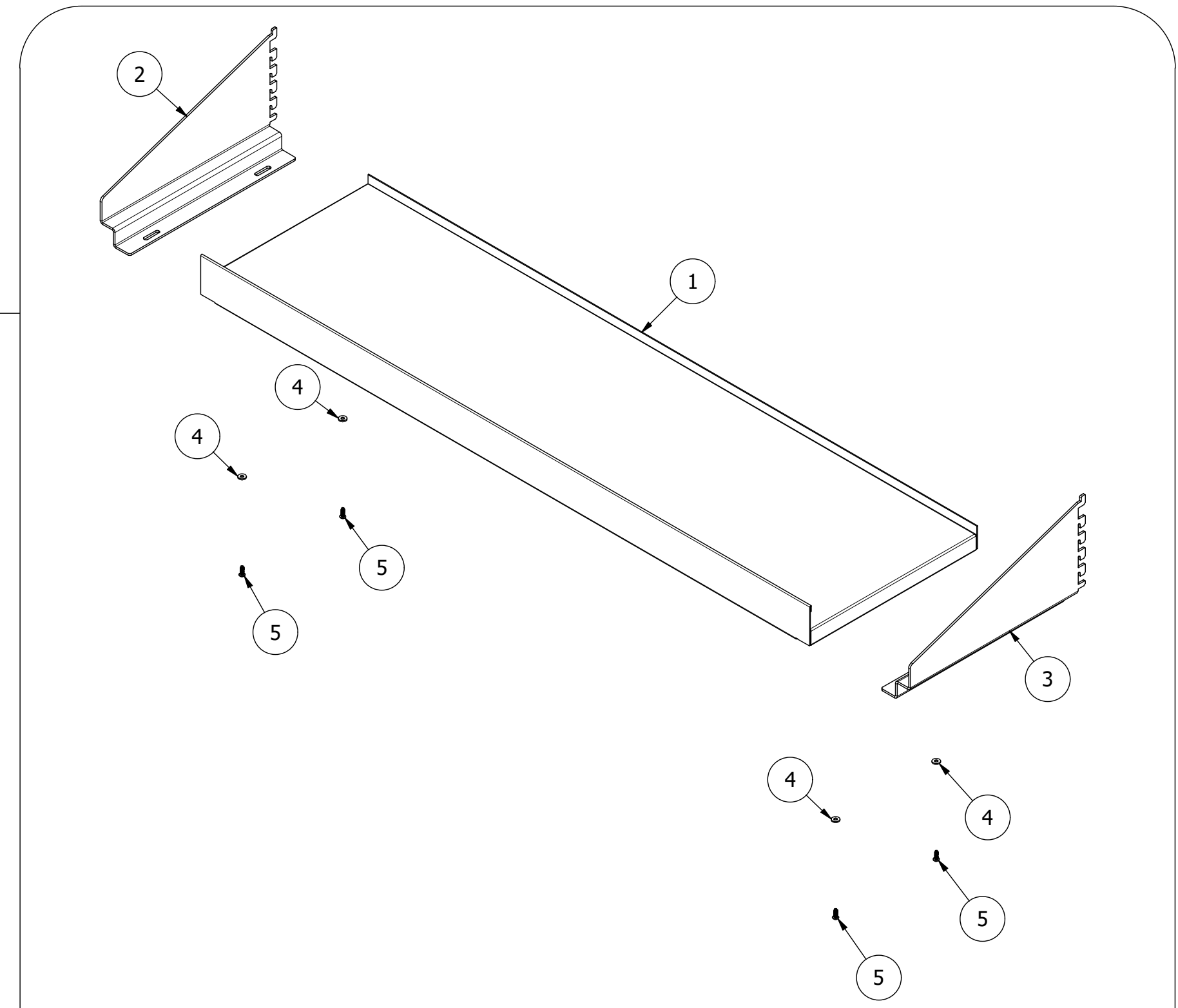
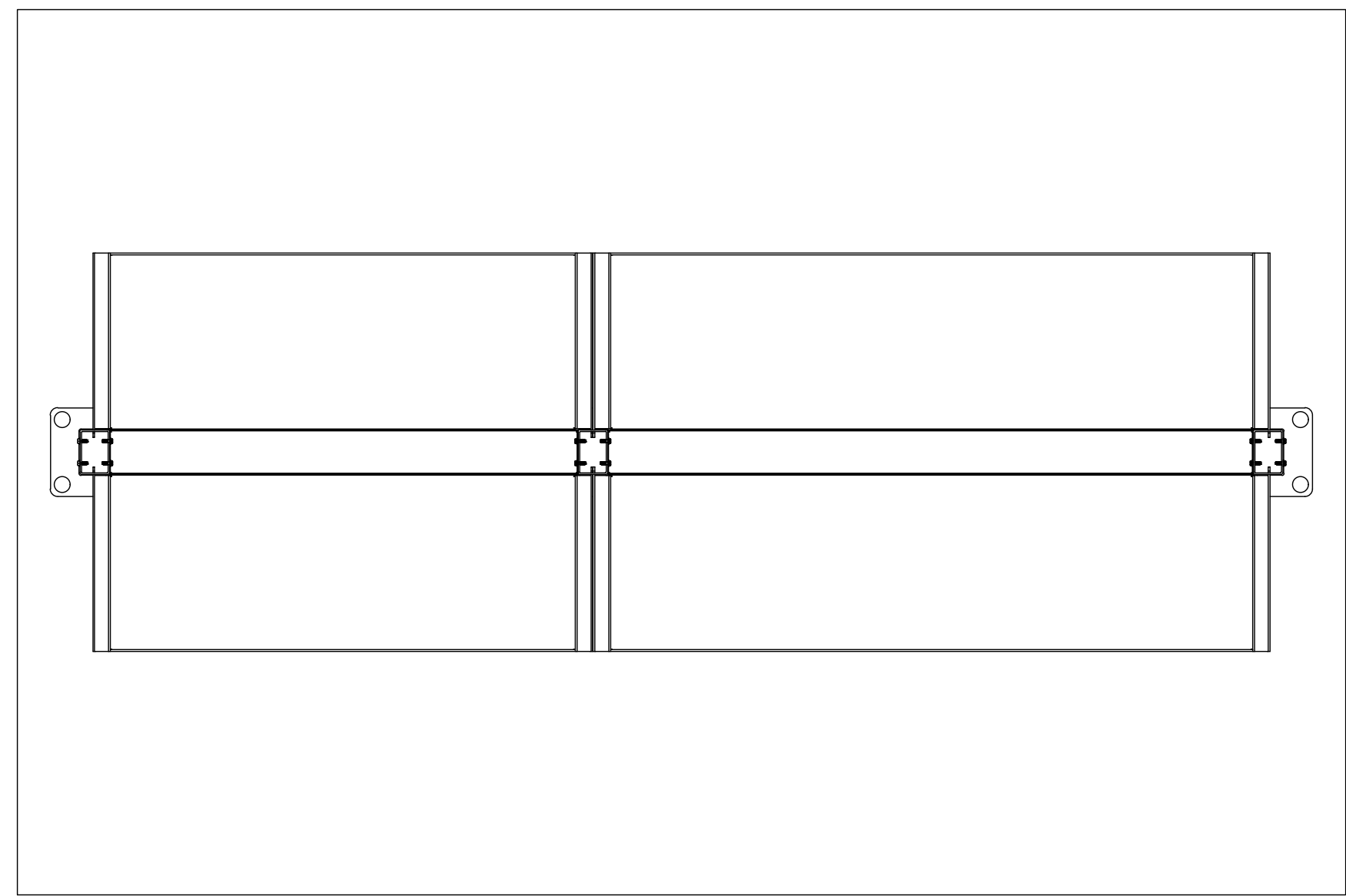


K13070-12 SHELF PARTS LIST			
ITEM	QTY	PART NUMBER	DESCRIPTION
1	1	C13070-12	SHELF ASSY, 31.5x12, G2
2	1	C13026-1	BRKT, SHELF, 12 LH, G2
3	1	C13026-2	BRKT, SHELF, 12 RH, G2
4	4	H236	SAE WASHER, #8
5	4	H138	SCREW, 8-32 THRD-CUT

K13032-12 PARTS LIST			
ITEM	QTY	PART NUMBER	DESCRIPTION
1	1	C13032-12	SHELF ASSY, 43.5x12, G2
2	1	C13026-1	BRKT, SHELF, 12 LH, G2
3	1	C13026-2	BRKT, SHELF, 12 RH, G2
4	4	H236	SAE WASHER, #8
5	4	H183	SCREW, 8-32 THRD-CUT

K13033-12 SHELF PARTS LIST			
ITEM	QTY	PART NUMBER	DESCRIPTION
1	1	C13033-12	SHELF ASSY 55.5X12, G2
2	1	C13026-1	BRKT, SHELF, 12 LH, G2
3	1	C13026-2	BRKT, SHELF, 12 RH, G2
4	4	H236	SAE WASHER, #8
5	4	H183	SCREW, 8-32 THRD-CUT

K13034-12 SHELF PARTS LIST			
ITEM	QTY	PART NUMBER	DESCRIPTION
1	1	C13034-12	SHELF ASSY, 67.5X12, G2
2	1	C13026-1	BRKT, SHELF, 12 LH, G2
3	1	C13026-2	BRKT, SHELF, 12 RH, G2
4	4	H236	SAE WASHER, #8
5	4	H183	SCREW, 8-32 THRD-CUT

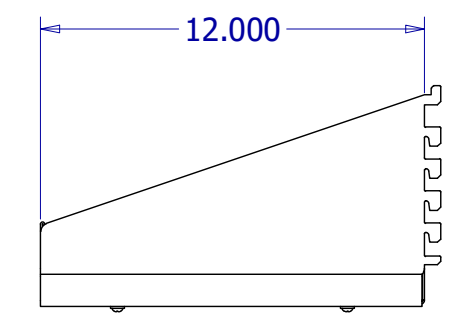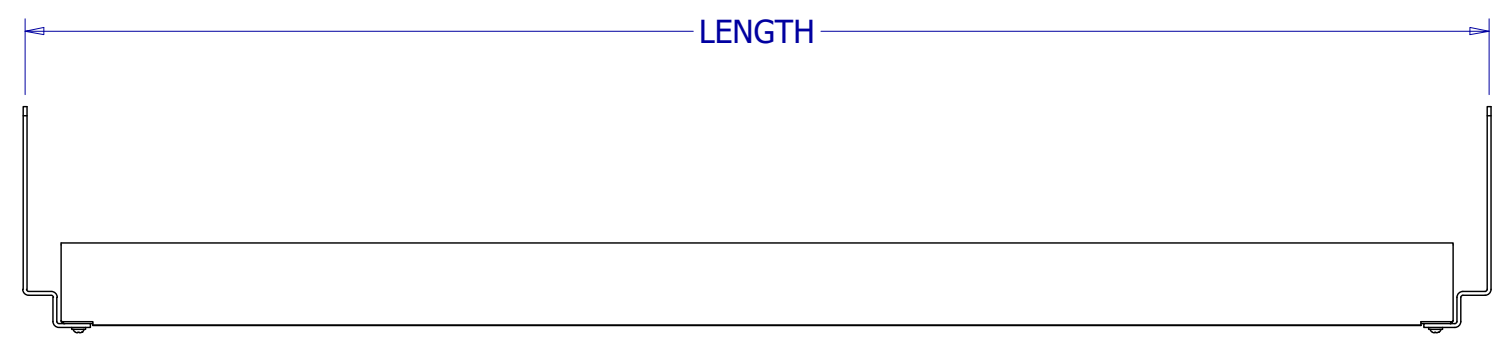

PART #	LENGTH
K13070-12	33.75"
K13034-12	45.75"



DRAWN AG	10/28/2022	LOOPED LOGIC 747 Calle Plano, Camarillo CA 93012 805-498-3121	
CHECKED		TITLE BENCHTOP UPRIGHT POST SYSTEM	
QA		FINISH PETAL WHITE	
APPROVED		SIZE D	PART NO
UNLESS OTHERWISE NOTED, GEN. TOLERANCE +/- .015"; ALL DIMENSIONS ARE IN INCHES	JOB # QUOTE #	SCALE 1 / 10	REV A
		MATL	SHEET 1 OF 1