

USER MANUAL

M-Series Laboratory Bench

M1, M2, M3, M4, M5, M6 & M-ERA



Version 1.4

HANSON LAB SOLUTIONS

747 Calle Plano, Camarillo CA 93012

(805) 498-3121 www.hansonlab.com



Content

<u>SAFETY INSTRUCTIONS</u>	02
<u>GROUNDING INSTRUCTIONS</u>	03
<u>CLEANING INSTRUCTIONS</u>	03
<u>ASSEMBLY INSTRUCTIONS</u>	03

Important Safety Instructions

When using this product, basic precautions should always be followed, including the following:

DANGER: TO REDUCE THE RISK OF ELECTRIC SHOCK:

- Always unplug this furnishing from the electrical outlet before cleaning.

WARNING: TO REDUCE THE RISK OF BURNS, FIRE, ELECTRIC SHOCK, OR INJURY TO PERSONS:

- Unplug from the outlet before putting on or taking off parts.
- Close supervision is necessary when these products are used by, or near, children, invalids, or disabled persons.
- Use these products only for their intended use as described in these instructions. Do not use attachments not recommended by the manufacturer.
- Never operate these products if they have damaged cords or plugs, if they do not work properly, or if they have been dropped or damaged. Return the products to a service center for examination and repair.
- Keep the cord away from heated surfaces.
- Never drop or insert any object into any opening.
- Do not use outdoors.
- Do not operate where aerosol (spray) products are being used or where oxygen is being administered.
- Max Load Rating for Countertops - 338 lbs.
- Max Load Rating for Upper Shelves – 169 lbs.
- Max Load Rating for Middle and Lower Shelves – 236 lbs.
- For loading always put heavier items at the bottom and not near the top to help prevent the possibility of the furnishing tipping over.

ELECTRICAL RATINGS: 120V, 60HZ, 20A PER CIRCUIT

CAUTION: Risk of Electric Shock - This unit has multiple power supply cords. Unplug all cords before moving or servicing this furnishing.

WARNING: Risk of Electric Shock - Connect this furnishing to a properly grounded outlet only. See Grounding Instructions.

SAVE THESE INSTRUCTIONS

Grounding Instructions

This product is rated more than 15 amperes and is for use on a circuit having a nominal rating of 120 volts and is factory-equipped with specific electric plug to permit connection to a proper electric circuit. Make sure that the product is connected to an outlet having the same configuration as the plug. No adapter should be used with this product.

This product may have more than one power supply cord. Ensure that all cords are unplugged before moving, testing or repairing the product.

Cleaning Instructions:

The table and light fixture can be cleaned with a damp cloth and mild cleaning solution. Solvents should not be used. If spray cleaner is used, spray on the cloth, not the furniture to avoid cleaning solution entering enclosed electrical areas.

Assembly Instructions:

Please refer to specific model number complete parts list and guide. Shelves will need to be installed in the following manner.

Assembling Shelves and Brackets: Shelves are attached to brackets with supplied screws using available holes.

Installing Shelves: Tilt the shelf and insert the upper hook in the upright slots first. While applying upward pressure, rotate the shelf until the lower hooks are fully inserted. Tap the shelf downward with a rubber mallet or a clenched fist. Ensure all hooks are fully engaged in the slots on the table upright.




For more information on the M-Series Bench Systems or for questions, please contact us.



 747 CALLE PLANO
CAMARILLO, CA 93012

 INFO@HANSONLAB.COM

 WWW.HANSONLAB.COM

 (805) 498-3121